



Laysan Bio Inc.

PRODUCT STORAGE INSTRUCTIONS

Because PEG reagents are temperature, air and moisture sensitive they should be stored in a freezer at or below -10C upon receipt. The Acrylate products are also light sensitive as well. Please allow PEG reagents to equal room temperature before opening. We recommend storing our PEG reagents as the dry, powder form only and not in any type of solution. Each time the bottle is opened it will need to be backfilled with an inert gas (i.e. Argon, Nitrogen) and tape the lid before being stored again. ACRL products should be stored in an amber or foil bag to prevent light exposure.

WE DO NOT RECOMMEND YOU STORE PEG PRODUCTS IN SOLUTION.

If your reagent requires rebottling please follow these steps:

Allow reagent to equilibrate to room temperature

While reagent is thawing prepare glove box or another suitable area for packaging which will allow for an inert atmosphere

Flush the repackaging enclosure with an inert gas (i.e. Argon, Nitrogen)

Repackage into desired aliquots

Flush each aliquot with inert gas before closing then tape lid and place in amber or foil bag to prevent light exposure

Store each aliquot at or below -10C until ready to use

LIGHT SENSITIVE

KEEP EXPOSURE TO LIGHT TO A MINIMUM

If you do not have the proper equipment to rebottle Laysan Bio offer a bottling service at \$50 set up fee and \$5.00 a bottle. This service must be ordered before the product ships.

If you have questions regarding this information we encourage you to contact our sales team at 256.586.9004 or sales@laysanbio.com.

For technical assistance email tech@laysanbio.com.

Laysan Bio Inc.
1560 Tower Drive
Arab, AL 35016
Phone: 256.586.9004
sales@laysanbio.com

1560 Tower Drive Arab, Alabama 35016

Phone: 256.586.9004 Fax: 256.586.9007